



MDF-U53VA-PA

VIP® Series

-86°C Ultra-Low Temperature Upright Freezer 18.3 cu.ft. | 519 L



Panasonic's VIP Series ultra-low temperature freezers are designed for safe, reliable storage of biologicals. Energy efficiency in VIP Series freezers is achieved through balanced temperature performance and power management. The advanced combination of cabinet design, electronics, refrigeration and critical components are engineered for enhanced security, performance and reliability.

Reliability You Can Count On

The robust refrigeration system is powered by Panasonic engineered compressors. The system provides dependable cooling by balancing temperature performance and energy management. Reserve refrigeration power assures exceptional temperature recovery following door openings. VIP freezers have demonstrated >99% reliability with excellent performance history.

Maximum Storage, Minimal Footprint

The upright freezer employs a patented VIP insulation composite that maximizes space efficiency and holds up to 352 2" boxes or 224 3" boxes within a 7.22 ft² (0.67 m²) footprint.

Uniform Preservation

The combination of patented VIP vacuum insulation panels with Panasonic designed compressors, specific to ultra-low temperature applications, assures superior uniformity (+/- 5°C) throughout the cabinet to maintain the integrity of the samples.



Microprocessor Controls/Alarms

Integrated microprocessor-based controller and LED display include comprehensive setpoint, alarms, monitoring, diagnostic and communication features.



Optimum Sample Storage

This space saving cabinet design offers a reduced wall thickness for maximum interior volume. Ideal for storage of DNA, RNA and other biological samples.



Extra Protection, Easy Access

The stainless steel framed, insulated inner doors with gasket seal firmly against the cabinet to help minimize change in interior temperatures during routine door openings. Inner door latches feature an ergonomic design for one-handed operation.

VIP -86°C ULT Upright Freezer

MDF-U53VA-PA




VIP Vacuum Insulated Panels

Inside-out engineering starts with a patented VIP insulation composite which results in an efficient, thin-wall cabinet. Insulated inner doors improve temperature uniformity, minimize cold air loss during door openings and extend warm-up times during power outages.

Model MDF-U53VA-PA Temperature Uniformity Data


-80°C setpoint at 20°C ambient temperature, empty freezer

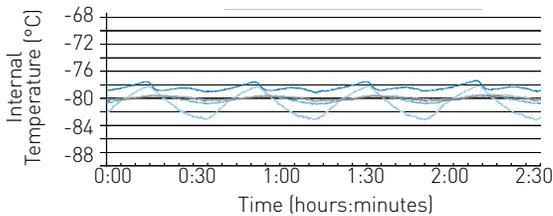

As interior air temperature stabilizes at the desired setpoint, mass storage volume sustains tight temperature uniformity of the load inside the insulated cabinet.

Dimensions

Unit : inch | mm

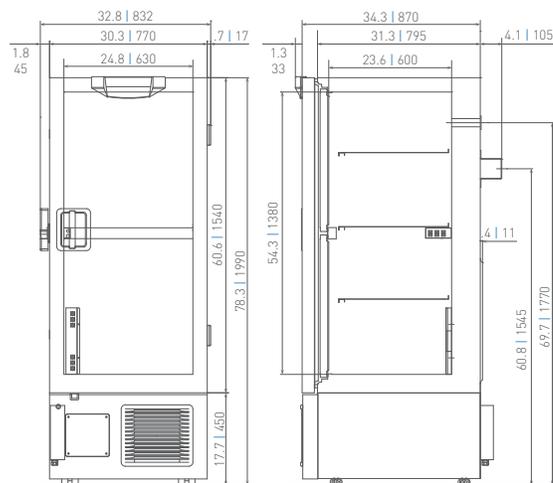

Model Number		MDF-U53VA-PA
External Dimensions (W x D x H) ¹⁾	inches mm	30.3 x 34.3 x 78.3 770 x 870 x 1990
Internal Dimensions (W x D x H)	inches mm	24.8 x 23.6 x 54.3 630 x 600 x 1380
Volume	cu.ft. liters	18.3 519
Net Weight	lbs. kg	639 290
Capacity (2" boxes)	qty	352
Capacity (3" boxes)	qty	224
Inventory Racks (2" boxes)	qty	16
Inventory Racks (3" boxes)	qty	16
Performance		
Warranty		Panasonic Five Year Warranty ²⁾
Cooling Performance ³⁾	°C	-86
Temperature Setting Range	°C	-50 to -90 in 1°C increments
Temperature Control Range ³⁾	°C	-50 to -86 in 1°C increments
Control		
Controller		Microprocessor, touchpad data entry with display lockout
Display		LED
Temperature Sensor		Pt-1000
Refrigeration		
Refrigeration System		Cascade
Compressors	W	(1) 450 High-Stage and (1) 750 Low-Stage
Refrigerant		CFC-free R-404A and R-508B
Insulation Material		Rigid polyurethane foam (PUF) + VIP
Construction		
Exterior & Interior Material		Painted steel
Outer Door	qty	1
Outer Door Lock		Standard
Inner Doors	qty	2 insulated, ABS, w/ stainless steel frame and positive latches
Shelves	qty	3
Shelf Dimensions (W x D)	inches mm	23.9 x 21.0 608 x 535
Max. Load - per Shelf	lbs kg	110 50
Max. Load - Total Freezer	lbs kg	440 200
Vacuum Release Port		Standard
Access Ports	qty	3
Access Port Position		Center upper back wall, left bottom and right bottom
Access Port Diameter	inches mm	0.6 17
Casters	qty	4 (2 leveling feet on front base)
Alarms		(V = Visual Alarm, B = Buzzer Alarm, R = Remote Alarm)
Power Failure		V-B-R
High Temperature		V-B-R
Low Temperature		V-B-R
Filter		V-B
Door Open		V-B
Electrical and Noise Level		
Power Supply		115V, 1Ø, NEMA 5-20P, requires NEMA 5-20R receptacle
Frequency	Hz	60
Noise Level ⁴⁾	dB (A)	51
Options		
Liquid CO ₂ Back-up System		CKV-UB21-PW
Temperature Recorder		
Circular Type	6", 7 day circular	MTR-C954 ⁵⁾
Chart Paper	52 charts per box	C7100386REV
Ink Pen	pack of 6	R252
Optional Communication System		
Wireless, cloud-based, automatic data management		LabAlert [®] Monitoring System



Panasonic Healthcare Corporation of North America
 1300 Michael Drive, Suite A, Wood Dale, IL 60191
 Toll Free USA (800) 858-8442, Fax (630) 238-0074
www.panasonic-healthcare.com/us/biomedical

¹⁾ Exterior dimensions of main cabinet only, excluding handle and other external projections

²⁾ Current warranty offered at time of printing and may be subject to change

³⁾ Air temperature measured at freezer center, ambient temperature +30°C, no load

⁴⁾ Nominal value - Background noise 20dB(A)

⁵⁾ Requires sensor cover MTR-DU700SF-PW